

Guest Medical

ALKALI NEUTRALISER POWDER



500g

ORDER CODE: H9424

Instructions for use:

USE WITH THE GUEST MEDICAL CHEMICAL SPILLS KIT

- Always wear personal protective equipment provided in the kit
- If possible, dilute caustic spillages with water before applying the Alkali Neutraliser Powder
- Apply powder to alkali spillage slowly from the outside to the centre
- The mixture will turn **yellow** when the **spillage is completely neutralised**
- If permitted, the mixture can be flushed away with water without damage to the environment
- Otherwise, collect granule/spillage mixture using disposable Guest Medical Hazard Scoops and Scrapers and discard into waste bag provided according to your organisational policy

Safety precautions:

- Avoid contact with Hydrofluoric acid
- Do not use on body fluid spillages

Manufacturer

Guest Medical Limited
Aylesford ME20 6SW, United Kingdom
T: +44 (0) 1622 791895
E: info@guest-medical.co.uk
W: www.guest-medical.co.uk

This quantity will neutralise;
500ml Sodium hydroxide 50%
600ml Potassium hydroxide 45%
1000ml Ammonium hydroxide 30%

CITRIC ACID, EC-NUMBER 201-069-1



WARNING

- Causes serious eye irritation
- Causes skin irritation
- May be harmful if swallowed
- Do not get in eyes, on skin, or on clothing
- Avoid breathing dust
- Wear protective gloves/eye protection

STORAGE

Store in original container.
Keep container tightly closed in a cool, dry, well ventilated place.

Batch No:

Expiry: